## ACURA® capillar 846

### POSITIVE DISPLACEMENT MICROPIPETTE DIREKTVERDRÄNGUNGS-MIKROPIPETTE MICROPIPETTE A DEPLACEMENT POSITIF

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#### SAFETY PRECAUTIONS

- Read operating instructions carefully; keep available for future reference.
- Observe manufacturer's instructions and guidelines for operation and maintenance.
- Refer to and follow regulations about handling of potentially hazardous reagents and glass components.
- Before use, check instrument for good working conditions.
- Do not use harsh chemicals (such as acetone) to clean instrument.
- This product should be used only for its intended purpose.
- Mind possible handfatigue during serial pipetting and its medical consequences (such as repetitive strain injuries RSJ).

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#### DESCRIPTION

The Acura® capillar 846 is a positive displacement micropipette line intended for accurate, precise and easy pipetting.

It has a wide range of applications such as:

- Viscous liquid distribution
- Dosing of volatile liquids
- Chemical and cosmetic applications
- Use of oils and detergents

Ajustable micropipette with 5 preset volumes. The plunger draws the sample into the precision glass capillary without any dead volume. No air interface means that volume dispensed will always be the same regardless of viscosity. The plunger is designed to deliver the sample without any noticeable residue, which allows the capillary to be reused with other fluids.

Five pipettes cover the full range of usual sample volumes from 1 to 200 µl.

Each pipette is supplied with a control quality certificate.

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#### DOSING INSTRUCTION

Volume selection is a matter of seconds. Simple mechanical device gives high precision, eliminate risks of errors. At all times the selected volume clearly appears on the barrel.

#### Volume setting (Fig. 1)

Barrel D will slide after rotating from locked position ▼ to selector position I. Place selected graduation mark (for example 50 µl) flush with end of ring C. Then lock into position ▼.

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## Dispensing

1. Depress plunger fully and immerse capillary into solution. Release slowly.
2. As needed, wipe carefully external of glass capillary without touching the opening.
3. Place capillary against wall of receiving vessel. Smoothly depress plunger until it stops. While moving the capillary along the vessel wall.

**Note:** For volumes of 1  $\mu$ l and 5  $\mu$ l, it is mandatory to eliminate any forming drop at end of capillary to insure good precision.

4. After work is completed, discard capillary or rinse in distilled water or saline solution by operating plunger rapidly a few times.

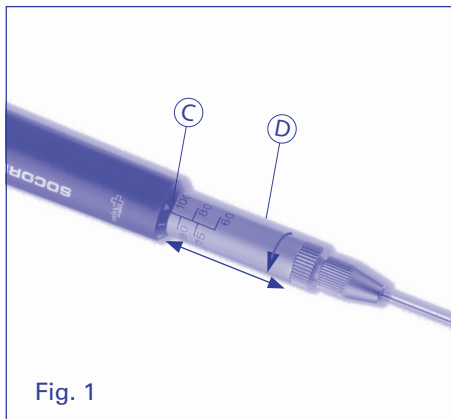
Fig. 1

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## MAINTENANCE

The instrument is designed to sustain long-term operation with minimal maintenance and service requirements.

Defective parts must be replaced, or instrument returned to dealer for repair. Order original spare parts from authorised dealers only.

The colour appearing on the finger rest of the pipette is to be found on the matching capillaries and spare plungers. Only assemble components with similar colour.

### Replacing and installing glass capillary (Fig. 2)

Loosen collet A and remove used capillary. Set pipette on **maximum volume** and insert new capillary **with colour mark first**.

**Do not try to remove a broken capillary over the plunger**, this would damage it. Remove plunger rod and broken glass over the rod. Remove collet A to eliminate any part of glass left in the collet A.

### Plunger replacement (Fig. 2)

To remove a worn plunger, loosen the collet A and carefully remove the glass capillary. Use the allen wrench to loosen the screw E. Remove the worn plunger.

Turn pipette upside down and insert a new plunger until it stops. Tighten screw E and reinstall capillary according to above.

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## CALIBRATION

Each positive displacement Acura® capillar 846 micropipette has been factory calibrated and controlled. The QC certificate accompanying the instrument includes control data and lot number. Calibration is carried out gravimetrically with bidistilled water at constant temperature ( $\pm 0.5^\circ\text{C}$ ) comprised between 20 and 25°C according to ISO 8655.

### Inlab calibration (Fig. 2)

Loosen collet A. Line up calibration mark B of capillary with the flat end of the plunger as shown in Fig. 2. Tighten collet A. The instrument is ready for operation. When plunger rod is fully depressed, the plunger tip will slightly protrude from glass capillary.

**Note:** Control with gravimetry method the desired precision for volumes < 50  $\mu\text{l}$

## WARRANTY

Your positive displacement Acura® capillar 846 micropipette is guaranteed against any material or manufacturing defects for the period of time specified in its QC certificate. Damages due to non-respect of manufacturer's instructions, safety precautions, glass breakage, as well as material colour alteration are excluded from the warranty. Repair and replacement of parts do not extend warranty time.

Should regular maintenance not eliminate a detected defect, return the instrument to the dealer from whom it was purchased after obtaining return authorization.

**Decontaminate instrument prior to returning it.**

## PERFORMANCE DATA AND ORDERING INFORMATION

Volumes $\mu\text{l}$	Inaccuracy (E%)			Imprecision (CV%)			Socorex Cat. No	Color Code
	Min. vol.	Mid. vol.	Max. vol.	Min. vol.	Mid. vol.	Max. vol.		
1-2-3-4-5 $\mu\text{l}$	< $\pm 3.0\%$	< $\pm 2.3\%$	< $\pm 2.0\%$	< $\pm 2.5\%$	< $\pm 2.0\%$	< $\pm 1.2\%$	846.005	White
5-10-15-20-25 $\mu\text{l}$	< $\pm 1.5\%$	< $\pm 1.2\%$	< $\pm 0.8\%$	< $\pm 1.0\%$	< $\pm 0.9\%$	< $\pm 0.6\%$	846.025	Grey
10-20-30-40-50 $\mu\text{l}$	< $\pm 1.2\%$	< $\pm 1.0\%$	< $\pm 0.7\%$	< $\pm 0.8\%$	< $\pm 0.7\%$	< $\pm 0.4\%$	846.050	Green
60-75-80-90-100 $\mu\text{l}$	< $\pm 0.7\%$	< $\pm 0.7\%$	< $\pm 0.7\%$	< $\pm 0.6\%$	< $\pm 0.5\%$	< $\pm 0.3\%$	846.100	Blue
100-120-150-175-200 $\mu\text{l}$	< $\pm 0.7\%$	< $\pm 0.7\%$	< $\pm 0.7\%$	< $\pm 0.3\%$	< $\pm 0.3\%$	< $\pm 0.3\%$	846.200	Red

\* Remove liquid meniscus from capillary end after aspiration, otherwise inaccuracy may be as high as 12%.

Performance values obtained with bidest. water at constant temperature ( $\pm 0.5$ ) comprised between 20 and 25°C in accordance with ISO 8655.

## ACCESSORIES

Spare glass capillaries Color code	Volume range	Packing	Socorex Cat. No.
White	1 – 5 $\mu\text{L}$	200 / pk	341.005
2 x white	5 – 25 $\mu\text{L}$	200 / pk	341.025
Green	10 – 50 $\mu\text{L}$	200 / pk	341.050
Blue	60 – 100 $\mu\text{L}$	200 / pk	341.100
Red	100 – 200 $\mu\text{L}$	100 / pk	341.200

Spare plungers with capillaries Color code	Volume range		
Steel (plunger alone)	1 – 5 $\mu\text{L}$	5 / pk	342.005
White	5 – 25 $\mu\text{L}$	5 / pk	342.025
Green	10 – 50 $\mu\text{L}$	5 / pk	342.050
Blue	60 – 100 $\mu\text{L}$	5 / pk	342.100
Red	100 – 200 $\mu\text{L}$	5 / pk	342.200

Work station for three pipettes Acura® capillar 846		1 / pk	320.340
---	--	--------	---------

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