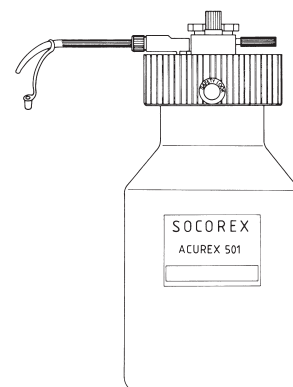
# ACUREX™ 501

## COMPACT DISPENSER KOMPAKTER DISPENSER DISPENSER COMPACT

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#### SAFETY PRECAUTIONS

- Read operating instructions carefully before use. Keep available for future reference
- Observe manufacturer's recommendations
- Follow regulations on handling of potentially hazardous reagents
- Before each use, check instrument for perfect working condition, in particular if glass parts have been broken.
- Before dosing, remove tip cap from delivery jet
- Prime and rinse dispenser carefully when left unused for a while or before disassembling
- Delivery jet should never point towards a person while dosing
- Never try to separate blocked parts by using force
- Cover delivery jet with cap, lock red safety screw and hold reagent reservoir firmly when carrying dispenser
- Maximal ambient temperature of use: < 50°C
- Do not use dispenser for purposes other than intended for
- Use only original spare parts

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#### DESCRIPTION

The Acurex™ 501 is a sturdy, uncomplicated precision dispenser with adjustable volume. All parts in contact with the liquid are of glass or FEP (no metallic springs) allowing safe handling of all reagents as well as diluted acids (except HF) and most solvents. The dispensing module is located inside the wide mouth amber glass reservoir for full protection. A retractable column gives lowest profile for convenient storage in a fridge.

Easy setting and fast priming account for good accuracy. Reproducibility is better than 0.5%. The unit can be fully disassembled for cleaning, all parts being interchangeable. It is autoclavable at 121°C fully assembled.

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## OPERATION

### Inspection and Installation

Examine unit for possible breakage; any defect should be reported to the dealer through whom the unit was purchased.

Carefully screw the delivery jet No. 13 – supplied unmounted – into the ETFE union located on the cap (see Fig. 1).

**Note:** Column stop No. 1 and upper stop No. 21 must be tightened before operating instrument.

### Safety

The dispenser head cannot be removed unless safety lock No. 30 has been loosened. For your safety, make sure "Safety Lock" is tightened after placing dispenser head on reservoir.

### Priming

Fill reservoir bottle with the desired reagent and place dispenser head on reservoir; gently rotate cap to obtain a tight fit. Loosen thumb screw No. 18 and slide indicator column No. 3 to largest volume; tighten thumb screw. Operate plunger handle No. 2 once on the full scale length and follow with several very short strokes to completely eliminate last air bubbles. The priming reagent may be returned to the reservoir if it has been dispensed into a clean vessel.

### Volume setting

When setting the volume, make sure that plunger handle No. 2 is pressed completely down. Slide column until desired volume appears at indicator ring No. 2b moulded on handle. Tighten thumb screw No. 18.

### Dispensing

The unit is now ready to operate; best results are obtained by moving the plunger gently and regularly between the upper and lower stops. Your movements will be correct if no droplet forms at the delivery jet.

After use, place tip cap over delivery jet to prevent withdrawal of the fluid and possible drying out of the ball valves which would result in deposition of undesired particles on the valve seat, thus impairing sealing efficiency.

## MAINTENANCE

### Disassembling

Make sure barrel contains no liquid by depressing plunger down to lower stop. Unscrew safety lock No. 30 and remove dispensing head from the reservoir. Prime if necessary.

Plunger can be easily removed by loosening thumb screw No. 18 then rotating upper stop No. 1 by 180°.

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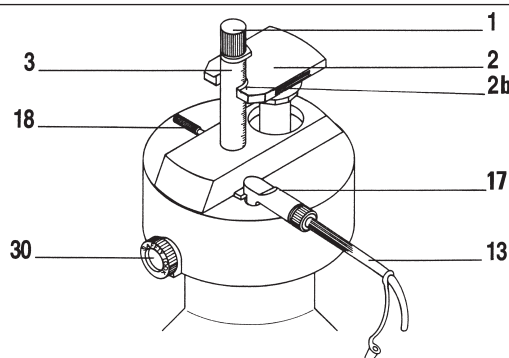
Fig. 1

## Cleaning

Clean fluid path regularly, i.e. whenever bottle has been emptied or before dosing another liquid.

When plunger is dragging or sticking, disassemble instrument. Soak dirty parts in a mild laboratory detergent and rinse carefully. To separate a frozen plunger, never use force, the glass barrel could break. Leave frozen plunger assembly soaking in an appropriate detergent overnight.

## Reassembling

Damaged parts should be exchanged, using only original spare parts. After reassembling, make sure all connections are tight and check for correct functioning.

## Calibration

No calibration required. Should performance no longer correspond to manufacturer's specifications, check for damaged parts or send to dealer for servicing.

## Sterilization

The Acurex™ 501 can be autoclaved at 121°C, 20 minutes (1 ATM) fully assembled, after priming and rinsing. Wait until instrument cools down before using it. Autoclaving also possible with PTFE extension tubing (cat. No. 1.523) Correct autoclaving and resulting sterility are the responsibility of the user.

## Materials

Parts in contact with liquid are chemically inert:

- Valve: Pyrex glass and synthetic ruby
- Barrel: neutral borosilicate glass
- Plunger: FEP coated glass
- Delivery jet and union: PTFE/ETFE/PFA
- External parts of dispenser: PPS/PP

**Note:** Screw No. 8 also available in PVDF (resistance to corrosive liquids)

**Warning:** Instrument not suitable for HF, concentrated strong acids and some specific solvents. Damages due to incompatible reagents are not warranty covered. In case of doubt, check for chemical compatibility or refer to manufacturer before use.

## PERFORMANCE DATA AND ORDERING INFORMATION

DISPENCER ACUREX 501/502					CAT. No. – KAT. Nr. – N° CAT.	
Volume Range Arbeitsvolumen Volume de travail	Division Teilung Division	Imprecision (CV%) Unpräzision (V%) Précision (CV%)			Reservoir Capacity Flascheninhalt Réservoir capacité	Reservoir with amber glass Flasche aus GW-Braunglas Réservoir verre jaune
		Min. vol.	Mid. vol	Max. Vol.		
0.2 - 2 mL	0.1	< 0.5%	< 0.35%	< 0.1%	250 mL	501.02025
0.2 - 2 mL	0.1	< 0.5%	< 0.35%	< 0.1%	1000 mL	501.021
0.2 - 2 mL	0.1	< 0.5%	< 0.35%	< 0.1%	2000 mL	501.022
0.4 - 5 mL	0.2	< 0.5% <sup>1)</sup>	< 0.35%	< 0.1%	500 mL	501.0505
0.4 - 5 mL	0.2	< 0.5% <sup>1)</sup>	< 0.35%	< 0.1%	1000 mL	501.051
0.4 - 5 mL	0.2	< 0.5% <sup>1)</sup>	< 0.35%	< 0.1%	2000 mL	501.052
1 - 10 mL	0.2	< 0.5%	< 0.35%	< 0.1%	1000 mL	501.101
1 - 10 mL	0.2	< 0.5%	< 0.35%	< 0.1%	2000 mL	501.102
1 - 30 mL	1.0	< 0.5% <sup>2)</sup>	< 0.35%	< 0.1%	2000 mL	501.302

<sup>1)</sup> at 0.5 mL    <sup>2)</sup> at 3 mL

Values obtained by a regular and smooth plunger movement with Aqua bidest. at constant temperature (±0.5°C) comprised between 20 and 25°C in accordance with ISO 8655.

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## WARRANTY

The Acurex™ 501 dispenser is guaranteed for a period of two years against any material or manufacturing defects. Damages due to non-respect of manufacturer's instructions, safety precautions or autoclaving conditions, as well as glass breakage are excluded from the warranty. Repairs and replacement of parts do not extend the warranty period.

Should regular maintenance not eliminate a detected defect, return the instrument to the dealer from whom it was purchased, after obtaining return authorization.

**Rince and decontaminate instrument prior to returning it.**

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