


Quick Guide

Profillette™ 406

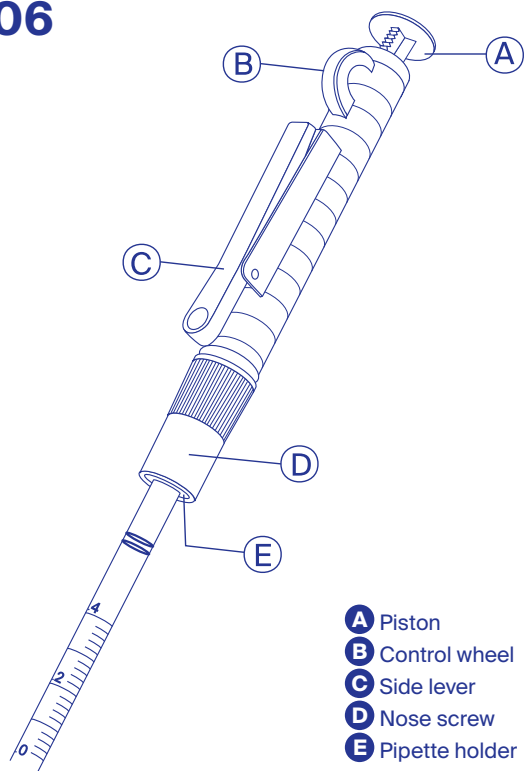

Related accessories	Packaging	Cat. No.
Spare silicone pipette holder	1 / pk	1.406.01
Replacement nose screw	1 / pk	1.406.02




Package contents

Pipette helper, Quick Guide

⚠ Safety precautions


- Read carefully and observe manufacturer's instructions and guidelines.
- Refer to and follow regulations about handling of potentially hazardous reagents.
- Before use, check instrument for good working conditions.
- Do not use harsh chemicals (such as acetone) to clean instrument.
- Do not use with liquids incompatible with silicone, PP, PTFE and ABS.
- Do not allow liquids to enter the instrument.
- Instrument is intended for liquid transferring in laboratory setting only.

Step-by-step instructions

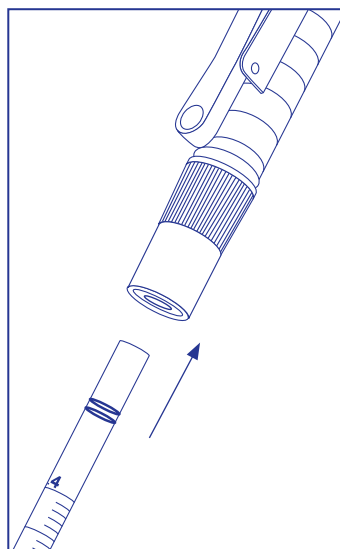


Pipette fitting

Carefully insert glass or plastic pipette into pipette holder (E).

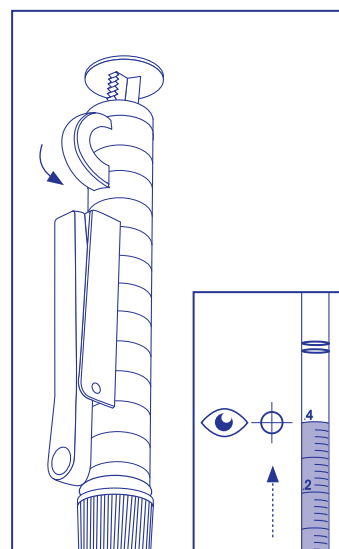


Aspiration

Hold vertically and rotate control wheel (B) to aspirate. Stop at slightly higher level than required volume, then adjust precisely.

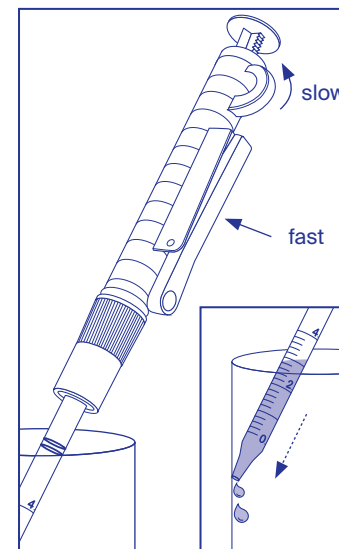

Dispensing

Hold pipette against recipient's wall and turn wheel in opposite direction to dispense drop by drop. To release liquid rapidly or to expel remainder liquid, press side lever (C).

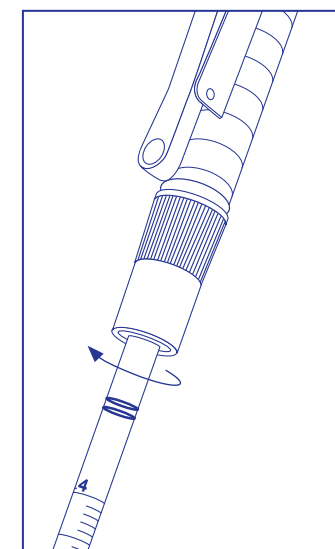

Pipette removing

Hold pipette and rotate gently pulling pipette from pipette holder.

Most common troubleshooting

Observation	Possible cause	Solution
Aspiration capacity is impaired	Silicone pipette holder damaged	Replace damaged parts
Pipette drips or pipette falls off	Pipette not properly positioned in silicone pipette holder	Verify pipette mounting
	Pipette holder or pipette damaged	Replace damaged parts

Scan to access operating instructions



EN / DE / FR